

Plastic Laminate Cabinetry

User's Guide



Mott Manufacturing LLC
562 Industrial Park Road
Maxwelton, West Virginia
USA, 24957
www.mott.ca



SAVE THESE INSTRUCTIONS

Table of Contents

Cleaning Pg.3

Notes of Caution Pg.3

Recommended Household Cleaners..... Pg.4

Back Panel Removal Pg.5

Drawer Removal Pg.6

Cleaning

To clean the surface, use a damp cloth or sponge and a mild soap. Rinse thoroughly with warm water and then wipe surface dry.

Difficult stains such as coffee can be removed by using a mild household cleaner and a soft bristled brush, repeat as necessary.

If a stain persists, use a baking soda and water paste. Apply and lightly scrub, excessive scrubbing or exerting too much down force could damage the laminate. So, proceed with caution.

Notes of Caution

1. Acidic or abrasive cleaners will more than likely damage laminate surfaces; do not use them.
2. Drain cleaners containing lye will permanently damage laminate surfaces. If you spill a drain cleaner, wipe it up immediately and rinse several times with water.
3. Hair, textile and food dyes can cause permanent stains. If dye should happen to spill, wipe it up immediately with an all-purpose cleaner.
4. Never place pots or pans directly from an oven or burner on an unprotected laminate surface; extreme heat could possibly cause blistering or cracking.

5. Do not work with oven cleaners, rust removers, or toilet cleaners near unprotected laminate. These items will damage the laminate. If a spill does occur, wipe up immediately and wash surface with soapy water and rinse several times.
6. Abrasive pads and steel wool will damage laminate. Do not use them for cleaning and do not store the steel wool on the surface of laminate, the metal will leave rust stains.
7. Prolonged exposure to bleach will cause discoloration
8. When cleaning laminate, DO NOT use products that contain the following ingredients:
 - Hydrochloric Acid
 - Sulfuric Acid
 - Hydrofluoric Acid
 - Phosphoric Acid
 - Sodium Hydroxide
 - Pumice

Recommended Household Cleaners

- Windex
- Glass Plus
- Fantastik
- Mr. Clean
- Formula 409

Removable Back Panels

To gain access to services behind cabinet:

- Remove shelf from cabinet
- Remove Philips head screws and pull out panel
- To replace back panel return to original position in cabinet and reinstall screws.



Drawer Removal

Full extension type drawer

- Pull the drawer out, all the way, push down on plastic tabs and pull drawer out to remove.
- To replace drawer, push full extension slide all the way into the cabinet.
- Line up channel on drawer housing with full extension channel in cabinet and push in. (Take care to line up runner channels as forcing them together may cause ball bearings to be dislodged and slide failure.)