

# TRI-ACCESS DOORS

# THE PIVOT™ SYSTEM



## Fully Accessible Door

The Pivot™ brings a new level of operator efficiency to floor mounted fume hoods with horizontal sliding sashes. Previously, all horizontal sliding sashes would only allow a 50% maximum opening for set up and tear down operation.

The Tri-Access Door System allows conversion from horizontal sliders to a hinged pivot that provides a 100% access opening; all done in 20 seconds with no tools! It's a major increase in efficiency offers a much quicker and safer access for users.



### Step 1

Disengage pin to open the hinge.



### Step 2

Fully slide the sash door to the hinged side before opening access door. Disengage the quick release latch and the door pivots to a fully open condition. Repeat with the second horizontal sash.



### Step 3

Lock sash with the bracket provided. Now larger items can be wheeled in at floor level.

**mott**  
GROUP OF COMPANIES

Steel Division & HQ - 452 Hardy Rd, Brantford ON, Canada N3T 5L8 • T. 519.752.7825  
Panel Products Division - 562 Industrial Park Rd, Maxwelton WV, USA 24957 • T. 304.497.2115

[www.mott.ca](http://www.mott.ca)

QUALITY BY DESIGN

2022.07